

RNAscope® VS Reagent Kit - BROWN

Description

The RNAscope® VS Reagent Kit – BROWN (Cat. No. 320600) is designed to work with the RNAscope® VS Target Probes (Cat. No. 300036) for fully automated *in situ* detection of RNA in formalin-fixed paraffin-embedded (FFPE) tissue specimens on slides using the Ventana™ DISCOVERY™ XT and ULTRA instruments. Each kit contains sufficient reagents in a convenient, Ready-To-Use (RTU) format to perform 60 slide/tests. The results of the assay can be viewed using a standard bright field microscope.

Contents and Storage

The components of the RNAscope® VS Reagent Kit – BROWN and quantities supplied are listed in the following table. The contents have a shelf life until the date of expiry indicated on the reagent label. It is important to follow the product storage conditions as specified on the label so that the reagents maintain their reliability.

Cat. No.	Sub-Kit	Component	Quantity	Storage
320043	RNAscope® Pretreat 2 Reagents	RNAscope® Pretreat 2	70 mL x 4 bottles	Room temperature (15–30 °C)
320060	RNAscope® VS Reagent Kit	RNAscope® VS Pretreat A	14 mL x 1 bottle	2-8 °C
		RNAscope® VS Pretreat B	14 mL x 1 bottle	2-8 °C
		RNAscope® VS AMP 1	14 mL x 1 bottle	2-8 °C
		RNAscope® VS AMP 2	14 mL x 1 bottle	2-8 °C
		RNAscope® VS AMP 3	14 mL x 1 bottle	2-8 °C
		RNAscope® VS AMP 4	14 mL x 1 bottle	2-8 °C
		RNAscope® VS AMP 5	14 mL x 1 bottle	2-8 °C
		RNAscope® VS AMP 6	14 mL x 1 bottle	2-8 °C
320630	RNAscope® VS Accessory Kit (complimentary kit)	RNAscope® VS Hematoxylin	7 mL x 1 bottle	2-8 °C
		RNAscope® VS Blueing Reagent	7 mL x 1 bottle	2-8 °C

IMPORTANT! Before running the assay, please download and review online the RNAscope® VS Reagent Kit-BROWN User Manuals (Cat. No. UM 320491 and UM 320493) at <http://www.acdbio.com/technical-support/downloads>.

Precautions

Handle biological specimens and other materials according to local laboratory guidelines and regulations. Avoid cross-contamination of samples and reagents to ensure quality results. Use equipment and reagents as specified by the manufacturer.

For safety and biohazard guidelines, refer to the “Safety” section in the RNAscope® VS Reagent Kit - BROWN User Manuals. For every chemical, read the Material Safety Data Sheets (MSDSs) and follow the handling instructions available at <http://www.acdbio.com/support/technical-doc/category/msds>. Wear appropriate protective eyewear, clothing, and gloves.

Contact Information

Address: 3960 Point Eden Way, Hayward, CA 94545 USA
Phone: 1-877-576-3636 option 1 **Fax:** 1-510-576-8798
Web: www.acdbio.com **Email:** support@acdbio.com

Copyright

©2015, Advanced Cell Diagnostics, Inc. All rights reserved.

Trademarks

RNAscope® is a trademark of Advanced Cell Diagnostics, Inc. VENTANA and DISCOVERY are trademarks of Roche. All other trademarks belong to their respective owners. Advanced Cell Diagnostics, Inc. reserves the right to change its products and services at any time to incorporate technological developments. This insert is subject to change without notice. Although this insert has been prepared with every precaution to ensure accuracy, Advanced Cell Diagnostics, Inc. assumes no liability for any errors, omissions, or for any damages resulting from the use of this information.