

RNAscope® Fluorescent Multiplex Reagent Kit

Description

Use the RNAscope® Fluorescent Multiplex Reagent Kit (Cat. No. 320851) along with the RNAscope® Probe – New Target Probe (Cat. No. 300031) for simultaneous *in situ* detection of up to three RNA targets in fresh frozen (FF) slide-mounted samples. Each kit contains sufficient reagents in a convenient, Ready-To-Use (RTU) format to perform 20 slide/tests. The results of the assay can be viewed using a fluorescent microscope with the appropriate filters.

Contents and Storage

The components of the RNAscope® Fluorescent Multiplex Reagent Kit and quantities supplied are listed in the following table. The contents have a shelf life until the date of expiry indicated on the reagent label. It is important to follow the product storage conditions as specified on the label so that the reagents maintain their reliability.

Catalog No.	Sub-Kit	Component	Quantity	Storage
320842	Tissue Pretreatment	RNAscope® Pretreat 4 (For Fresh, Fixed Frozen Tissues)	4 mL x 2 bottles	2-8 °C
		RNAscope® Pretreat 3 (cultured cells, PBMCs)	4 mL x 1 bottle	2-8 °C
320851	RNAscope® Fluorescent Multiplex Reagents	RNAscope® AMP 1-FL	3 mL x 1 bottle	2-8 °C
		RNAscope® AMP 2-FL	4.5 mL x 1 bottle	2-8 °C
		RNAscope® AMP 3-FL	3 mL x 1 bottle	2-8 °C
		RNAscope® AMP 4-FL-Alt A Display module	4.5 mL x 1 bottle	2-8 °C
		RNAscope® AMP 4-FL-Alt B Display module	4.5 mL x 1 bottle	2-8 °C
		RNAscope® AMP 4-FL-Alt C Display module	4.5 mL x 1 bottle	2-8 °C
		RNAscope® DAPI	3 mL x 1 bottle	2-8 °C
310091	RNAscope® Wash Buffer Reagents	RNAscope® Wash Buffer (50x)	60 mL x 4 bottles	15-30 °C

IMPORTANT! RNAscope® Reagent Kits share the same RNAscope® Wash Buffer, but have unique RNAscope® Detection Reagents and RNAscope® Pretreat Reagents based on different tissue types. Do not interchange the reagent components of the Detection Reagents, even those having the same name. Refer to the latest versions of user manuals Part 1 for Sample Preparation and Part 2 for the RNAscope® Detection Assay available online at <http://www.acdbio.com/technical-support/downloads>.

Precautions

Handle biological specimens and other materials according to local laboratory guidelines and regulations. Avoid cross-contamination of samples and reagents to ensure quality results. Use equipment and reagents as specified by the manufacturer.

For safety and biohazard guidelines, refer to the "Safety" section in the *RNAscope® Fluorescent Multiplex Reagent Kit User Manual* (Cat. No. UM 320293). For every chemical, read the Material Safety Data Sheets (MSDSs) and follow the handling instructions available at <http://www.acdbio.com/technical-support/downloads/technical-doc/category/msds>. Wear appropriate protective eyewear, clothing, and gloves.

Contact Information

Address: 3960 Point Eden Way, Hayward, CA 94545 USA
Phone: 1-877-576-3636 option 1 **Fax:** 1-510-576-8798
Web: www.acdbio.com **Email:** support@acdbio.com

Copyright

©2015, Advanced Cell Diagnostics, Inc. All rights reserved.

Trademarks

RNAscope® is a trademark of Advanced Cell Diagnostics, Inc. All other trademarks belong to their respective owners. Advanced Cell Diagnostics, Inc. reserves the right to change its products and services at any time to incorporate technological developments. This insert is subject to change without notice. Although this insert has been prepared with every precaution to ensure accuracy, Advanced Cell Diagnostics, Inc. assumes no liability for any errors, omissions, or for any damages resulting from the use of this information.